


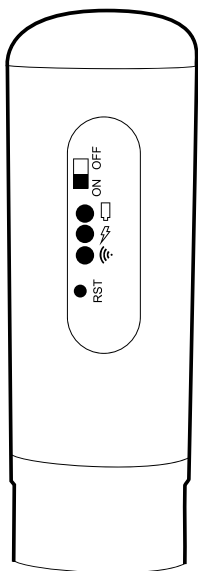

MICRO i

Digital Wi-Fi Imager
For IOS & Android Mobile Devices




Parts Of The Imager	3
Set Up	4
Using The Application	6
Internal Battery	6
Reset Button	6
Technical Data	7
Warranty	8

Notes Before Use

- Do not disassemble the product.
- Do not clean the product with alcohol organic solvents.
- Do not touch the lens with your fingers.
- Avoid outdoor use.
- If the product gets wet, leave PC connection immediately. Do not disassemble or try to dry product. Send to repair center if the product was affected by liquid or other elements.
- Measurement data is only for reference.
- Fully charge product before using. When charging, the red LED light will be off automatically indicating that the internal battery is fully charged.
- Do NOT use product while charging.
- Portable power supply can also be used to charge the product.
- It is normal for the product to get hot with continuous use of 10 minutes. Turn microscope off and allow it to cool off for 30 mins before re-using.

Parts Of The Imager

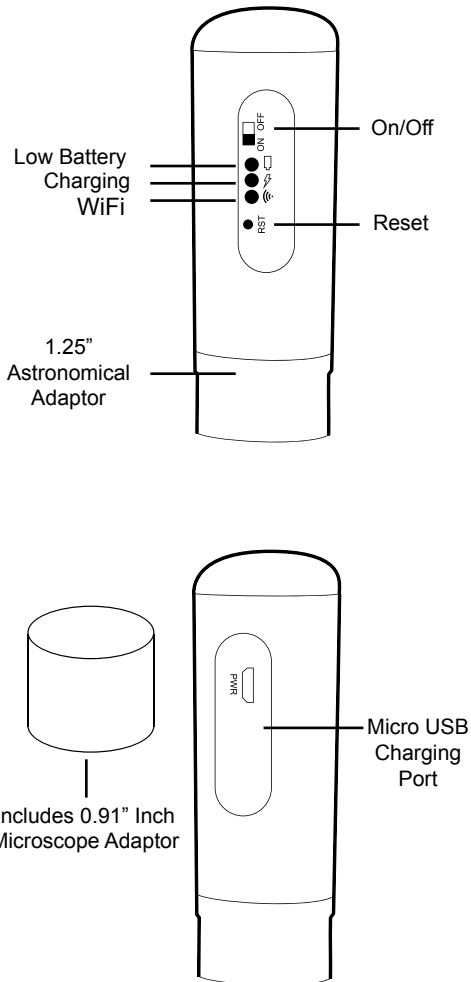

Set Up

Note For Android Users

You will need to change your settings to allow installation of apps from unknown sources



1. Turn on the microscope

Scan QR code to download and install the WiFi Microscope application to run on device.

Andriod Devices



IOS Devices





Once installed pair the devices

For cellphone

1. Go to Settings.
2. WiFi search for BARSKA Microscope Imager and select and link to BARSKA-Microscope-Imager_305038.
3. Enter Password: 12345678 to connect the devices.
4. Select the WiFi scope application to run the program.

Using The Application

Place appropriate adaptor into eyepiece.

 <p>Take Images</p>	<p>Press the camera button and save file.</p>
 <p>View Files</p>	<p>Press the folder button. Choose image or video option to view file.</p>

Internal Battery

Product can NOT be used when the red LED Low Battery Indicator light is on.

Connect product via USB port to a power source. The red LED Low Battery Indicator will be off once the product is fully charged.

Reset Button

If program malfunctions, use a small tool to press the reset button to restart product.

Technical Data

<p>Focus Range</p>	<p>0 mm - 40mm Other Ranges Can Be Customized WiFi Standard : IEEE 802.11 b/g/n Standards Compliant</p>
<p>Image</p>	<p>Image 640*480 CMOS Sensor</p>
<p>Software</p>	<p>WiFi Capable IOS and Android</p>
<p>Application Software</p>	<p>Wireless Microscope.apk/ APP Illumination Scope: 0 - 30000LUX</p>
<p>Language</p>	<p>WiFi Capable IOS and Android - English, other languages can be customized</p>
<p>Power</p>	<p>Rechargeable 1500mA Self-Protection Lithium Battery Charging Time 3 Hours Continuous Use 2.5 Hours</p>
<p>Specs</p>	<p>Height: 150 mm Width: 40 mm Weight: 110 g</p>
<p>Storage Temperature</p>	<p>0°C - 40°C</p>
<p>Humidity</p>	<p>45% RH - 85 RH%</p>
<p>Range</p>	<p>49 feet</p>

1 YEAR LIMITED WARRANTY

MICROSCOPE

BARSKA warrants your microscope to be free from defects in materials and workmanship for one (1) year. BARSKA will repair or replace such product or part thereof which, upon inspection by BARSKA, is found to be defective in materials or workmanship. As a condition to the obligation of BARSKA to repair or replace such product, the product must be returned to BARSKA together with proof-of-purchase satisfactory to BARSKA.

The proper Return Merchandise Authorization Number (RMA) must be obtained from BARSKA in advance of return. Please e-mail info@barska.com or call 1.888.666.6769 for Return Merchandise Authorization number (RMA) before any returns. NOTE: All merchandise received without a valid RMA number will be returned to shipper at his/her own expense.

BARSKA products should be sent to the following address for warranty repairs. Products must be packed carefully and sturdily to prevent damage in transit, and returned freight prepaid to:

BARSKA
855 Towne Center Drive
Pomona, CA 91767

For additional and updated information please visit our website at www.barska.com

Please include all of the following when returning BARSKA products for service and/or replacement:

1. Please write your complete details (Name, Address, Telephone#, E-mail address, RMA#, etc.)
 2. Purchase receipt or Proof of Purchase. (Original/Copy)
 3. A Brief explanation of the defect.
 4. A Check/Money Order for \$25.00 to cover inspection, shipping and handling.
- *Please allow 6-8 weeks for delivery.

The customer shall be responsible for all costs of transportation and insurance, both to and from BARSKA, and shall be required to prepay such costs.

BARSKA shall use reasonable efforts to repair or replace any microscope covered by this warranty within 30 days of receipt. In the event repair or replacement shall require more than 30 days, BARSKA shall notify the customer accordingly. BARSKA reserves the right to replace any product which has been discontinued from its product line with a new product of comparable value and function.

This warranty shall be void and of no force of effect in the event a covered product has been modified in design or function, or subjected to abuse, misuse, mishandling or unauthorized repair. Further, product malfunction or deterioration due to normal wear is not covered by this warranty.

BARSKA DISCLAIMS ANY WARRANTIES, EXPRESS OR IMPLIED, WHETHER OF MERCHANTABILITY OF FITNESS FOR A PARTICULAR USE, EXCEPT AS EXPRESSLY SET FORTH HEREIN. THE SOLE OBLIGATION OF BARSKA UNDER THIS LIMITED WARRANTY SHALL BE TO REPAIR OR REPLACE THE COVERED PRODUCT, IN ACCORDANCE WITH THE TERMS SET FORTH HEREIN. BARSKA EXPRESSLY DISCLAIMS ANY LOST PROFITS, GENERAL, SPECIAL, INDIRECT OR CONSEQUENTIAL DAMAGES WHICH MAY



RESULT FROM BREACH OF ANY WARRANTY, OR ARISING OUT OF THE USE OR INABILITY TO USE ANY BARSKA PRODUCT. ANY WARRANTIES WHICH ARE IMPLIED AND WHICH CANNOT BE DISCLAIMED SHALL BE LIMITED IN DURATION TO A TERM OF ONE YEAR FROM THE DATE OF ORIGINAL RETAIL PURCHASE.

Some states do not allow the exclusion or limitation of incidental or consequential damages or limitation on how long an implied warranty lasts, so the above limitations and exclusions may not apply to you.

This warranty gives you specific legal rights, and you may also have other rights which vary from state to state.

BARSKA reserves the right to modify or discontinue, without prior notice to you, any model or style microscope.

If warranty problems arise, or if you need assistance in using your microscope contact:

BARSKA Customer Service Department Tel. 1.888.666.6769 Fax. 909.445.8169 Monday-Friday 8:30AM-5:00PM PST or e-mail: service@barska.com


NOTE: This warranty is valid to U.S.A. customers who have purchased this product from an authorized BARSKA dealer in the U.S.A. This warranty is limited to the original purchaser and is not transferable.

BARSKA shall not be liable for any consequential, incidental and/or contingent damages whatsoever and is not responsible for damage to personal property. We will not pay shipping, insurance or transportation charges from you to us, or any import fees, duties and or taxes. This warranty supersedes all previous BARSKA warranties.



MICRO i

Digital Wi-Fi Imager
For iOS & Android Mobile Devices

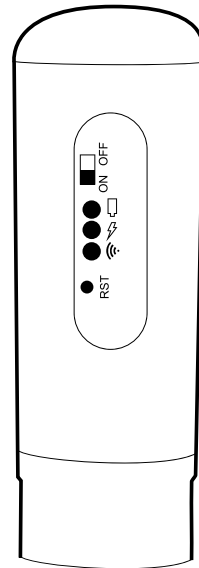

All trademarks and copyrights are the property of their respective owners. Featured brand trademarks are not associated with this product.